



Protein Standards for Mouse Angiogenesis ELISA Strip

Catalog Number EA-1022

(For Research Use Only)

Protein Standards


<u>Tube</u>	<u>Protein</u>	<u>Stock Conc.</u>	<u>Volume</u>
#1	Mouse TNFa	400ng/ml	10ul
#2	Mouse IGF-1	400ng/ml	10ul
#3	Mouse VEGF	400ng/ml	10ul
#4	Mouse IL-6	400ng/ml	10ul
#5	Mouse FGFb	400ng/ml	10ul
#6	Mouse IFNr	400ng/ml	10ul
#7	Mouse EGF	400ng/ml	10ul
#8	Mouse Leptin	400ng/ml	10ul

Preparation of protein standard dilutions

1. Add 200ul of Diluent buffer the wells of the first strip, and add 100ul of Diluent buffer to the wells of the rest strips according to the following table.
2. Add appropriate amount of diluted protein standards to the first strip according to the following table 1.
3. Use multi-channel pipette to mix the dilutions and transfer 100ul to the next dilution wells.
4. Repeat the transfer until 5th strip. Trash 100ul from 5th strip after mixing. Do not transfer to the 6th Strip

Note: Substrate incubation time may vary due to different antibodies reactivity. Stronger signals (Strong blue color) could be stopped early after 5 minutes. Weaker signals should be incubated for 10-30 minutes. Always stop protein standards along with samples from the same row at the same time.

Table 1: Dilution of protein standards

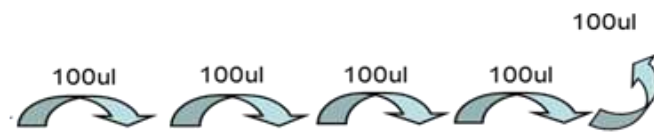

<u>Protein</u>	<u>Added to 1st Strip</u>	<u>1st Strip</u>	<u>2nd Strip</u>	<u>3rd Strip</u>	<u>4th Strip</u>	<u>5th Strip</u>	<u>6th Strip</u>
Mouse TNFa	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse IGF-1	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse VEGF	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse IL-6	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse FGFb	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse IFNr	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse EGF	4ul	200ul	100ul	100ul	100ul	100ul	100ul
Mouse Leptin	4ul	200ul	100ul	100ul	100ul	100ul	100ul
			1:2	1:4	1:8	1:16	Blank

Table 2: Concentrations of protein standard dilutions

<u>Protein</u>		<u>1:2</u>	<u>1:4</u>	<u>1:8</u>	<u>1:16</u>	<u>Blank</u>
Mouse TNFa	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse IGF-1	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse VEGF	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse IL-6	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse FGFb	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse IFNr	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse EGF	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank
Mouse Leptin	8ng/ml	4ng/ml	2ng/ml	1ng/ml	0.5ng/ml	Blank